

user manual

pc_o.1300



pc_o.
imaging

In this manual you find instructions for the pco.1300 camera in its standard version for the end-user and the modified pco.1300solar version.

Target Audience: This camera is designed for use by technicians, engineers, and scientists.

In case of any questions or comments, please contact us at PCO.



telephone	+49 (0) 9441 2005 50
fax	+49 (0) 9441 2005 20
email	info@pco.de
postal address	PCO AG Donaupark 11 93309 Kelheim, Germany

The cover photo shows an exemplary PCO camera system.
The lens is sold separately.

Copyright © 2010 PCO AG (called PCO in the following text), Kelheim, Germany. All rights reserved. PCO assumes no responsibility for errors or omissions in these materials. These materials are provided "as is" without warranty of any kind, either expressed or implied, including but not limited to, the implied warranties of merchantability, fitness for a particular purpose, or non-infringement. PCO further does not warrant the accuracy or completeness of the information, text, graphics, links or other items contained within these materials. PCO shall not be liable for any special, indirect, incidental, or consequential damages, including without limitation, lost revenues or lost profits, which may result from the use of these materials. The information is subject to change without notice and does not represent a commitment on the part of PCO in the future. PCO hereby authorizes you to copy documents for non-commercial use within your organization only. In consideration of this authorization, you agree that any copy of these documents, which you make, shall retain all copyright and other proprietary notices contained herein. Each individual document published by PCO may contain other proprietary notices and copyright information relating to that individual document. Nothing contained herein shall be construed as conferring by implication or otherwise any license or right under any patent or trademark of PCO or any third party. Except as expressly provided, above nothing contained herein shall be construed as conferring any license or right under any PCO copyright. Note that any product, process, or technology in this document may be the subject of other intellectual property rights reserved by PCO, and may not be licensed hereunder.

Updated August 2010 © PCO AG

Table of Contents

1	Introduction	4
	The advantages of the pco.1300 for your success	
2	Safety Instructions	5
	Save money, reduce stress, stay healthy	
3	System Components	6
	What's in the Box?	
4	Installation	7
	Get your camera and computer ready for start.	
5	Quick Start	8
	How to get an image fast and easy	
6	Camera Control for the pco.1300	9
	The main interface for all camera settings	
	6.1 The "Camera Control" window	9
	6.2 The <i>Timing</i> sheet	10
	6.3 The <i>Sensor (Size)</i> sheet	11
	6.4 The <i>Sensor (Misc.)</i> sheet	12
	6.5 The <i>CamRam</i> sheet.....	14
	6.6 The <i>Recording</i> sheet	14
7	Camware Features	15
	Software for image acquisition and archiving	
	7.1 If no camera is connected	15
	7.2 Overview	16
	7.3 The File Menu.....	17
	7.4 The Camera Menu.....	21
	7.5 The Acquisition Menu.....	22
	7.6 The View Menu.....	23
	7.7 The Window Menu.....	25
	7.8 The Help Menu.....	26
	7.9 The Local Menu.....	27
	7.10 Additional Features.....	28
Appendix		
A1	Technical Data	30
A2	Mechanical Dimensions	31
A3	Pinout of 15pin DSUB Connector	32
A4	Image File Formats	33
A5	Service and Maintenance	35
A6	Customer Service and Trouble Shooting	36
A7	Firmware Update	37
A8	Digital Camera Tools CD-ROM	38
	About PCO	39

1 Introduction

pco.1300: when you need a scientific grade camera with a compact design for system integration as well as end-user applications.

Features

This versatile high performance 12 bit CCD camera system has an extraordinary quantum efficiency of up to 65%. At the heart of the camera is an FPGA processor allowing for sophisticated control and accurate timing of the CCD and associated electronics. In addition, a proprietary offset control algorithm has been developed which provides very high offset stability, regardless of ambient temperature or signal changes, ensuring accurate and repeatable quantitative data over long periods of time.

The camera's main features are:

- excellent resolution (1392 x 1040 pixel)
- superior quantum efficiency up to 65%
- cooled 12 bit dynamic range
- superior low noise of $6e^-$ rms @ 10MHz
- no glow functionality at long exposure times
- excellent DSNU at long exposure times
- optimal offset stability and control (<1count)
- hot pixel correction integrated
- cooling of -25°C vs. ambient
- standard IEEE1394a interface

OEM


The pco.1300's most unique feature is its flexibility for customization to fit any OEM user application. ROI, binning, cooling, as well as custom housing, submicron alignment, and other features of the camera can be selected and optimized to accommodate the user's application.

End-user

In its standard version this camera serves as a well-priced alternative for the scientist, whose focus is on sophisticated low noise performance rather than on fair design. **This manual is about this standard version.**


pco.1300

This low light camera system is perfectly suited for many sensitive and low noise imaging applications, such as biochip reading, fluorescence imaging, confocal & light microscopy etc.

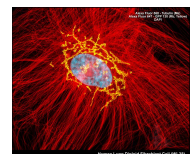

pco.1300solar

This modified standard version is particularly designed for electroluminescence (EL) applications in the range of 850nm to 1200nm.

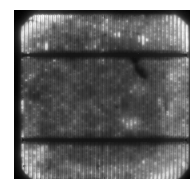

2 Safety Instructions



Never operate the camera in humid or dusty environments or in places with high amounts of X-ray radiation. **Humidity, dust or X-rays could damage the camera.**



To avoid the risk of water condensation, protect the camera against extreme changes of ambient temperature. **If condensation enters the camera, there is the risk of electric shock.**



To prevent damage to the camera, **the system must be kept stable and protected against strong jolts or vibrations.** The socket at the bottom of the camera is to be used for mounting purposes only.



The slits in the camera case (side & back planes) are designed for heat dissipation by the camera fan. **To prevent overheating of the camera, do not block these slits.** Do not leave the camera system in direct sunlight to avoid the risk of overheating.



Electric shock warning – **Never slide any items through the slits into the camera** because of the risk of electric shock if the voltage parts inside are touched.



Each time the camera is used, **check the power cable for any damage.**



Never position the cable in a way that it could become a **tripping hazard.**



Do not force the lens onto the camera. To protect the lens connector thread from damage, use minimal force when attaching a lens to the camera.

If any of the following conditions apply, immediately switch off the camera, separate it from the power line and contact our customer support:




- ⚠ If the power cable or the power plug seems to be worn or damaged.
- ⚠ If liquids have penetrated the device.
- ⚠ If, after thoroughly reviewing the instruction manual, the device is still not operating properly.
- ⚠ If the camera has been dropped or the casing is damaged.

3 System Components

The camera system includes the following parts.

Camera Head

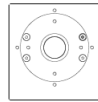

C-mount optical connection:

The distance between the front edge of the c-mount and the sensor is 17.52 mm. For standard c-mount lenses and adapters.



15pin DSUB connector for power and trigger input (connect to power supply)

IEEE1394a FireWire connector (connect to PC)

LEDs indicate camera status

The LED lights flash during initialization.

LED status	
red	camera busy
green	exposure



1/4-20 UNC and **3/8-16 UNC** thread for mounting



Serial Number tag



Trigger & Power Connection

Proprietary combined DSUB/BNC connector.

DSUB: for pinout see appendix A3

BNC: external exposure trigger in (TTL) (see chapter 6.2)



Power Supply

24V AC adapter with IEC14 chassis plug and proprietary DSUB/BNC connection cable



Power Cord

Standard IEC13 connector (for Germany only)



FireWire Cable (optional)

IEEE1394 FireWire cable 6pin/6pin (1.8m)



FireWire PC Card (optional)

32bit and 33MHz IEEE 1394 FireWire host adapter



Digital Camera Tools (CD-ROM)

The accompanying CD contains (see appendix A8):

- Camware: software for camera control & image acquisition
- camera driver & tools
- software development kit (SDK) and demo programs in C and C++

4 Installation

You will find all necessary files on the accompanying CD. You may also download the newest versions of our software, camera drivers, and third party software drivers from our website (support section).

Your pco.1300 is a special modification of a **pco.camera**.

Minimum system requirements:

- Clock speed > 1.6GHz
- RAM > 512MB
- Windows XP/ Service Pack 2 or Win2000/ Service Pack 4
- 1280 x 1024 pixel resolution display
- 32-bit-graphics card (16.7 million colors)

4.1 FireWire 400 (IEEE1394a)

Please insert the FireWire host adapter into the PCI expansion slot of your computer, if not already installed!

The drivers for FireWire interfaces or cards are part of Windows XP or part of the board manufacturer package. The installation of an additional hardware driver for the interface is not necessary.

4.2 Camera Driver

The first time the camera is connected via FireWire to the computer, Windows will notify the user that a new device has been detected. Afterwards, the typical Windows hardware drivers' installation process will begin. At this point, you are asked for the location of the corresponding *.inf file. Please click on the drive where the CD has been inserted or to the directory where the installation software has been downloaded. Windows will perform any further installation and system registration automatically.

4.3 Camware

The Camware 32-bit Windows application software enables you to control every camera parameter or setting. Images can be displayed on a monitor and may be downloaded and stored.

The CD runs automatically (if not: please begin the installation manually by double clicking on the setup.exe file). The installation software copies all necessary DLL files into the corresponding Windows directories. While doing this, older versions of these files will be replaced, if present. All required Windows registration entries will be made automatically. After a successful installation, you will find the program file 'Digital Camera Toolbox' in your program directory. Other helpful tools are also installed in the same file (see appendix A8).

To uninstall the Camware program, please use the *Software* feature under Window's *System Control*.

➤ **After installation, please restart/reboot your computer!**

5 Quick Start

Note: In order to get familiar with your new camera & software it might be helpful, if you first aim at an object that is easy to focus and that can be seen at standard light conditions.


5.1 Preparation




- ✓ installation is finished (see chapter 4)
- ✓ an appropriate lens is attached (remove cap!)
or the camera is attached properly to the microscope, spectrograph or other scientific device
- ✓ camera is connected to the PC


- ✓ computer is on
- ✓ camera is connected to the power supply and ready (LEDs off)

5.2 Start

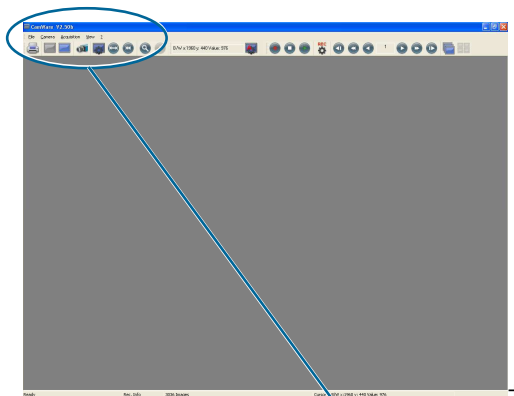- start Camware 



- open view window
- if not already open - 
- start "live preview" 
- apply "auto range peak" 
- you may adjust aperture and focus

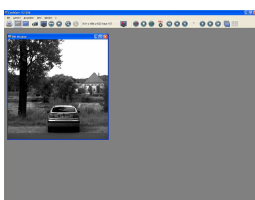

You should now clearly see the object in the window.

- If you need to change exposure time (e.g. the image is still either too dark or too bright), please go to chapter 6!
- If you want to record and save images, please see chapter 6 and chapter 7 for detailed information!

6 Camera Control for the pco.1300

The 'Camera Control' window in Camware is the main interface for all camera settings.


For further Camware features please see chapter 7!

6.1 The "Camera Control" window

The camera control window can be opened by selecting the proper command in the "Camera"-Menu or by the corresponding button in the toolbar:



The camera control dialog always adapts to the camera type connected. For the pco.1300 the camera control settings are spread over five **property sheets** ①, which will be explained below.

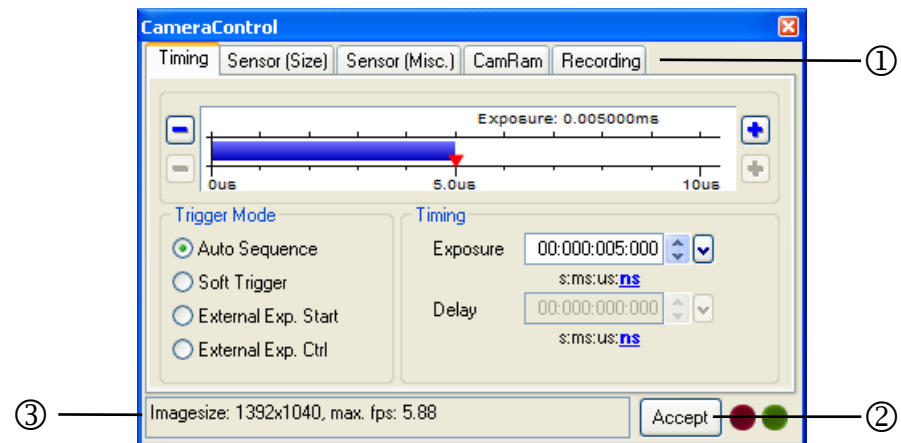

Changes to the camera control sheets must be completed by pressing the **"Accept" button** ②.

If the "Accept" button is not pressed, these changes will be ignored and lost. If the changes are within the allowed condition limits of the camera, a message is sent to the computer, the button disappears and the green lamp will be highlighted. If an illegal change was attempted, the camera will not accept the change and the red lamp will be highlighted.

In the **information field** ③ you can view some information about the camera. Click into the info field in order to scroll through the values.

pco.1300.solar.noglow.x2 Camera (1392x1040), b/w, no doubleshutter


El. Temp.: 25 °C, CCD Temp.: 9 °C, Power Temp. 0 °C

Imagesize: 1392x1040, max. fps: 10.52, est. rec. time: 00min 01s

Power Temp. not applicable for pco.1300.

Estimated recording time (est. rec. time) not applicable for pco.1300.

6.2 The “Timing” sheet

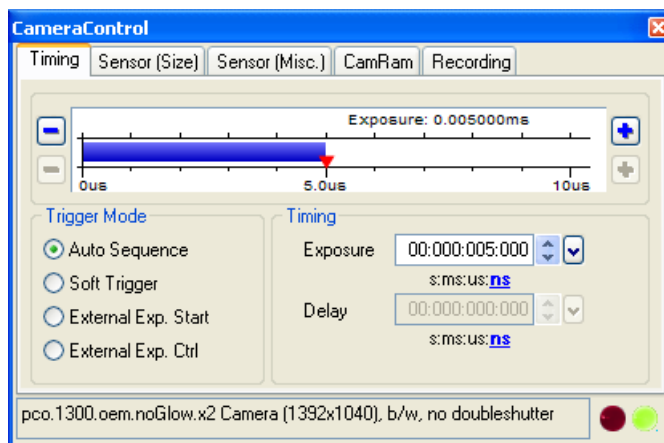

Trigger Mode

[Auto Sequence] The camera will optimize the image recording to achieve the best possible frame rate.

In the [auto sequence] exposure control mode, the camera determines the fastest possible image uptake rate depending on the adjusted exposure time and the required readout time. After a start command is given, the sequential recording is started until a stop command is given.

[Soft Trigger] Single images can be recorded with Camware control. The single image can be acquired by pressing the "Single Trigger" button. This button appears after pressing the "Start Record" button (see 7.2). Other signals cannot influence this operating mode.

[External Exp. Start] The image acquisition is triggered by an external signal. It is also possible to force a software trigger for a test image with the "Single Trigger" button.

In the [External Exp. Start] exposure control mode, single image recording is started by the falling or rising edge of the voltage signal at the BNC input (see chapter 3).

pixel clock	intrinsic + system delay
10 MHz	10 ... 160 µs
20 MHz	10 ... 100 µs

[External Exp. Ctrl] not available for pco.1300

Timing


The exposure time (delay time is not available for pco.1300) can be adjusted in steps of 158µs (10MHz) / 85µs (20MHz). For IR mode see page 12.

	normal mode	IR mode
pco.1300	5µs ... 1h	5ms ... 1h
pco.1300solar	5µs ... 1h	100ms ... 1h

In Camware exposure settings are possible in 5µs steps. However, the actual timing of the camera will be set to the nearest possible timing according to the 158µs (10MHz) / 85µs (20MHz) specification.

For example: 45µs setting in Camware will result in 5µs exposure, 50µs setting in Camware will result in 90µs exposure (@ 20MHz).

6.3 The “Sensor (Size)” sheet

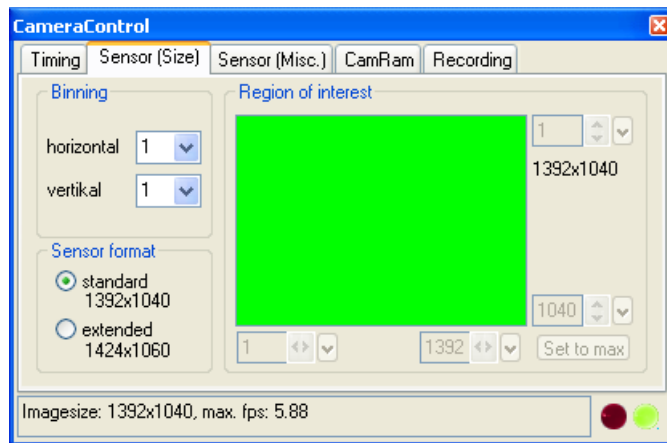

Binning

Binning combines neighboring pixels (in either the horizontal or vertical direction) to form super pixels. It increases the light signal of the resulting pixels and decreases the spatial resolution of the total image, which is recorded. Vertical binning (x2) will also double the speed.


Region of Interest

The ROI (region of interest) selects only a part of the sensor to be read out, in order to speed up the frame rate and to save memory (not available for the pco.1300).

Sensor Format

The implemented sensor has the option to readout a standard size (effective pixels recommended by the sensor manufacturer) or an extended size, which adds optical black pixels, which can be helpful for calibration or control purposes. In extended mode, you will not increase image size!

6.4 The “Sensor (Misc.)” sheet

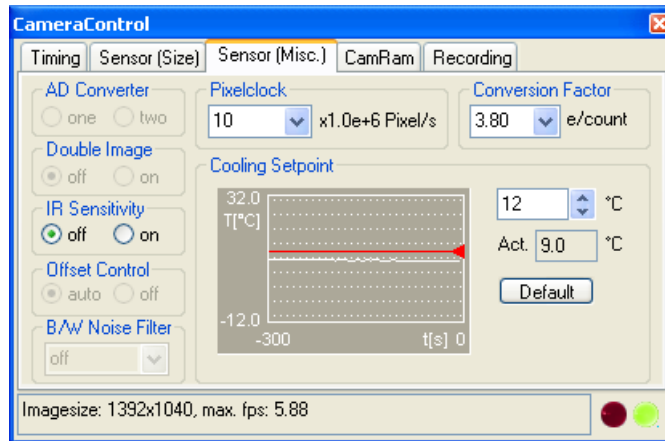

Pixelclock

The pixel clock sets the clock frequency and therefore the image sensor readout speed. At 10 MHz the image quality will be higher due to very low readout noise (read out at a lower bandwidth). At 20 MHz the image sensor is read out with double speed, achieving higher frame rates.

	10 MHz	20 MHz
pco.1300	5.9 fps	10.5 fps
pco.1300solar	5.9 fps	11.7 fps

AD Converter


Feature not available for pco.1300.

Double Image

This feature is widely used for particle image velocimetry (PIV) measurements, but not available for pco.1300.

IR Sensitivity = Low Light Mode

This feature uses a special image sensor control method, allowing greater sensitivity in the near infrared spectral range (red line in graph), in most cases at the expense of decreased anti-blooming performance.

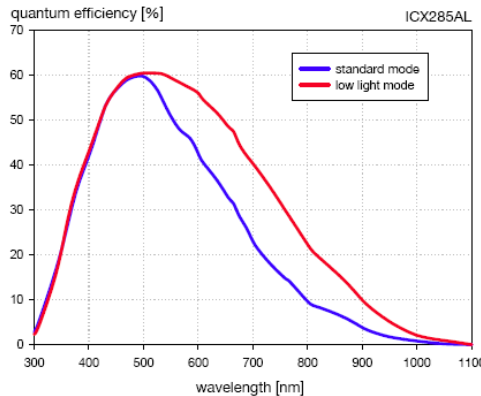

Offset Control

Offset control is done automatically for optimal results. User control is disabled.

B/W Noise Filter

In the pco.1300 an optimal hot pixel correction is already integrated. For exposure times > 1s you might need extra Hot Pixel List (see page 21).

Conversion Factor

Using the values from the dropdown list, the corresponding conversion factor can be selected. The conversion factor defines how many charges (electrons), which have been generated by light in the image sensor in each pixel, are necessary to generate one count (one intensity level) in the digital image. Therefore, the conversion factor describes the gain that is applied to the signal before it is converted into a digital value.


	selectable conversion factors	
pco.1300	3.5	3.9
pco.1300solar	1.9	3.8

Cooling Setpoint


The cooling set point adjusts the sensor temperature that should be reached by the camera system. The user should not set the cooling set point to the lowest possible temperature. If ambient temperature drifts, the controller will not have any overhead available for adjustment. The window shows the readout temperature of the image sensor and the adjusted temperature. The lower the temperature, the less dark current is accumulated, which is only important for longer exposure times. If exposure times are in the ms range or shorter, the lower temperature has no major influence. To keep the offset as constant as possible, moderate cooling is sufficient. When the "Default" button is pressed, it sets the cooling temperature back to the default setting.

We guarantee cooling of -25°C vs. ambient. In practice, cooling is possible of up to -40°C vs. ambient.

For the pco.1300solar only moderate cooling ($5^{\circ}\text{C} - 15^{\circ}\text{C}$) is preferable. Detailed quantum efficiency (QE) measurements revealed, that this is an advantage for EL applications. The graph clearly shows, that the QE increases for higher wavelengths at higher temperatures.

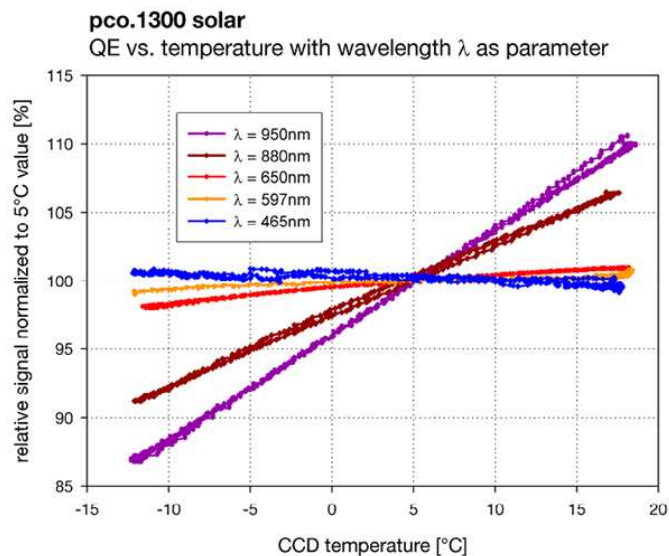


6.5 The “CamRam” sheet

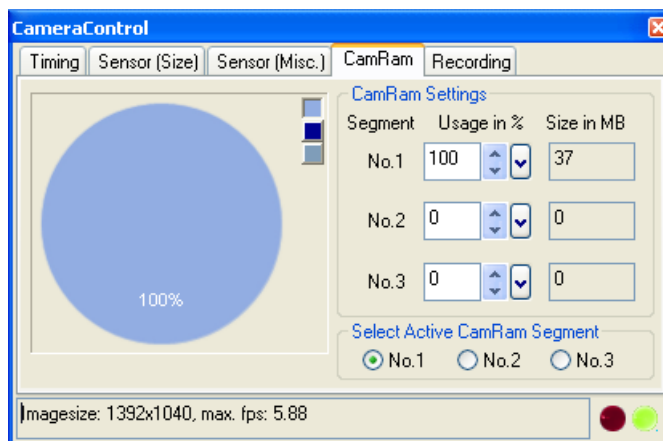

Feature not available in the pco.1300.

6.6 The “Recording” sheet

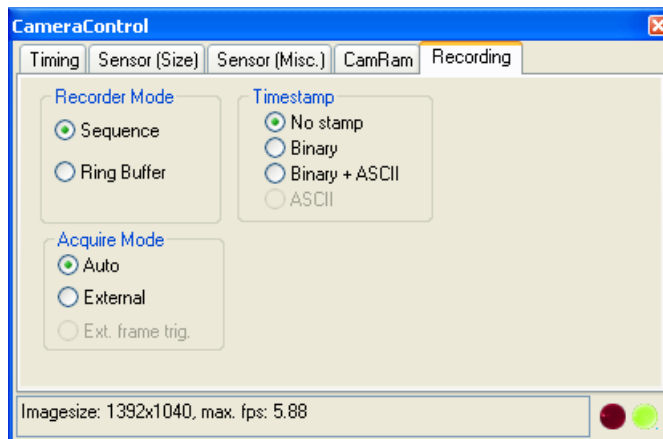

Feature not available in the pco.1300.

Recording with the pco.1300

The Camware will use free disk space on your computer. The recorded images will be temporarily saved as a 16bit multi TIFF under documents&settings\user\local\temp\PCO\cam0. Therefore, the camera runs in a virtual ‘Sequence’ Recorder Mode. The maximum number of images depends solely on the disk size of the computer.


Every new recording or shut-down of Camware will delete this temporary folder (including images)!

For further details on recording and saving images see the following chapters: 7.2; 7.3; 7.5!

Note: Multiple cameras can be connected with Camware, but only one pco.1300 camera can record! You may run two cameras simultaneously, but record with one after another.

7 Camware Features

PCO's renowned software for camera control, image acquisition, and archiving of images in various file formats.

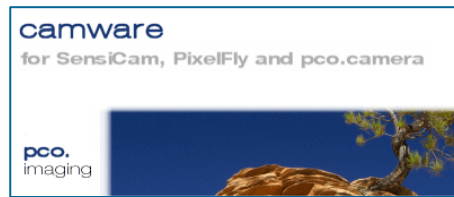


You will find an **online help** under the common Help menu.

Note: There are some features explained in the online help that are not applicable for this camera version.

7.1 If no camera is connected

When Camware is started, it automatically recognizes the camera type, which is connected to the computer and switched on.

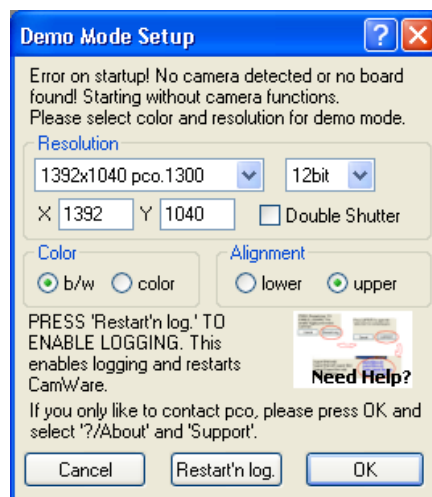

If Camware is started with no camera connected to the PC or with cameras switched off, it starts in demo mode.

In this mode all **image processing** features are available. The user only has to tell Camware which type of images will be opened. For that purpose, the "Demo Mode Setup" window opens and asks for the corresponding input.

Need Help? If this window pops up because you have trouble running the camera, please see instructions in appendix A6.

Resolution

The drop down list offers the existing image sensor spatial resolutions of all PCO camera systems. Please select the specific resolution and bit depth of the images to be opened! If double shutter images have been recorded and should be opened, "Double Shutter Mode" should be checked.

Color

With the radio buttons, the user can specify whether the image type is monochrome or color.

Alignment


These two radio buttons adjust whether MSB aligned ("upper") or LSB aligned ("lower") images have been stored (see page 18, "File Settings").

Settings to view the b16 files of the pco.1300:

	Resolution	bit	DoubleShutter	Color	Alignment
pco.1300	1392x1040	12	<no>	b/w	upper

7.2 Overview

Menu Bar

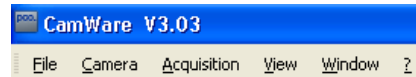

See chapter 7.3 - 7.8.

Local Menu



The “mouse right-click” is explained in chapter 7.9.

Toolbar



	prints the currently displayed image or comment window
	opens a b/w color display window
	opens the camera control window
	opens the convert control
	auto range peak/crop (see chapter 7.6 under “Convert Control BW”)
	live preview: useful for fast and easy camera adjustment and focusing This command disables the camera control window if it is opened!
	soft trigger: captures a single image (see chapter 6.2)

Cursor display




Col x:1640 y: 596 R:242 G:255 B:255	cursor: pixel position and intensity value
	continuous “auto range crop” scaling

Recorder



	record START / STOP / PLAY
	opens the recorder settings window
	replay navigation
 	The 'Recorder Browser' gives an overview of the recorded images. A click on the desired image makes it the current image in the Black/White or Color window. 'Skip #images' allows for convenient overview at large numbers of images.
	With the "Multi Frame Window" it is possible to view all active camera images in one window. The user might select between "time" oriented or "camera number" oriented order.

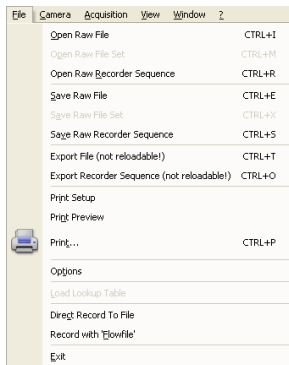

7.3 The File Menu

Open Raw File (single image only)

This command should be used to import a single image into the currently active image window. Only files with the extension and format of "*.b16" (PCO proprietary binary image format) and ".tif" (TIFF image format) can be imported. If the recorder is enabled, each imported image will be transferred to the buffer shown in the picture number. The image itself will be fitted to the current image size. If the recorder is disabled, the current image sizes will be set to the parameters of the imported image.

Open Raw File Set (for hsfc pro camera)

This command is used to open a set of images, previously saved with the 'Save Raw File Set' command. The corresponding windows will be opened automatically. Additionally a comment window will be opened, which shows the settings of the channel and a comment. This option is not accessible in case of the recorder is switched on.

Open Recorder Sequence (image sequence from one camera)

This command is used to import a sequence of images. If more than one camera is connected and an image window is currently open, the sequence will be loaded to the active window. If no image window is open, the images will be loaded to camera #1. This command opens the "Open file" dialog box. Only files with the extension and the format of "*.b16" and ".tif" (TIFF16) can be imported.

Save Raw File (single image only)

The "Save" command should be used to save the image, which is displayed in the active window. The command opens the "Save file" dialog box. The image file can be saved in 16bit "*.b16" and ".tif" format.

If more than one camera is connected, it is possible to save all current images by selecting "Export all images" in the "Save file" dialog box. With this feature it is possible to save one image of each active camera within one process step (it is not necessary to repeat the save process for each camera). The "Save" command will not be available, if there is no image window open. See 6.5 for further information!

Save Raw Set (for hsfc pro camera)

To save one image of each active camera within one single b16 file.

Save Raw Recorder (image sequence from one camera)

The "Save Recorder" command should be used to save or export image records. If more than one camera is connected and an image window is currently open, the record of the active window will be saved. The command opens the "Save file" dialog box. If another image format than "*.b16" or ".tif" (16bit) is chosen, image content will be lost, and it will be impossible to reload the images. The "Save Recorder" command will not be available, if there is no image window open. See 6.5 for further information!

Export File (not reloadable!)

Use this command to export the image in the active image window. This command will open the Save file dialog box. Files with the extensions "fts", "tif", "bmp", "asc", "jpg", and "jp2" can be exported. This item will be not visible, if there is no image window open.

Note: be aware of the different storage abilities of the formats, for example "*.bmp" - the bitmap format only allows for 8bit values to be stored and therefore the image content of a 12 or 14bit image is reduced, if stored as bitmap.

Export Recorder Sequence (not reloadable!)

Use this command to export a sequence of images. If more than one channel is connected and an image window is currently open the record of the window which has got the input focus will be saved. If no image window is open the images of channel 1 will be saved. This command will open the Export image dialog box. Files with the extensions "fts", "tif", "bmp", "asc", "avi", "mpg", "jpg", "jp2", and "wmv" can be exported.

If you export recorder files you will not be able to reload them into CamWare!

Print Setup

The "Print Setup" command opens the Windows "Print Setup" dialog box for adjustments and settings of the connected printer(s).

Print Preview

The "Print Preview" command opens a "Print Preview" window.

Print

Opens the Windows "Print" dialog box for the currently displayed image or comment display.

Options

This command opens the "Options" dialog, allowing special settings.

Options: File Settings



Tiff Bit Alignment

Here you can select the alignment of the pixel value. If this item is checked each pixel value is shifted from bit 12-1 to bit 16-5. This removes the 'dark image effect' inside other image applications.

Tiff RAW

Here you can select whether the 16bit RGB tif file is saved with 16bit raw or white balanced data. If this item is checked, each 16bit RGB pixel value is saved as raw 16bit RGB (48bit) value, without white balancing. Saving to 16bit RGB can be done by selecting the color display window, export recorder, selecting 8bit-tif (!) and checking 'Save 16bit RGB'.

ASCII file separator


Here you can select an appropriate separator for the ASCII file format.

Jpeg High Quality

Here you can select the quality of the saved jpeg/jp2 files. High quality saves images with 100% quality. Unchecking this option will create jpeg files with 85% quality.

Preserve Last Record

If checked, a message box will pop up in case the user may overwrite or loose data.

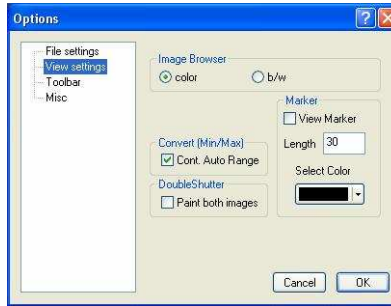Options: View SettingsImage Browser

Here you can select between showing all image tiles inside the browser window either b/w or color. Setting b/w will result in a faster refresh. (valid only for color CCD).

Marker


Use this command to display a cross in the centre of each image.

Convert (Min/Max) (“Continuous Auto Range Crop”, see chapter 7.6 under “Convert Control BW”)

This option enables the automatic min/max function during record and replay. Disable this feature if the application stops responding during this mode (due to operating system capacity problems).

Double Shutter:


This option enables the update of both double shutter images during record and replay. Disable this feature if the application stops responding during this mode (due to operating system capacity problems).

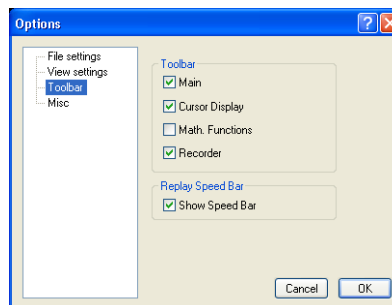
Options: ToolbarToolbar


Opens the corresponding toolbars (see 7.2)



Math. Functions

With the math window you have the ability to build the difference between a reference image and the actual image. If you activate math every new image acquired will be subtracted from the reference image or vice versa.

 last acquired image will be copied to reference buffer

 a reference picture will be acquired and copied to reference buffer


 add offset here to avoid negative values, which would not be visible

  enable / disable math function

Replay Speed Bar

The faster replay speed is useful if large sets of image data need to be scanned.

Options: Misc

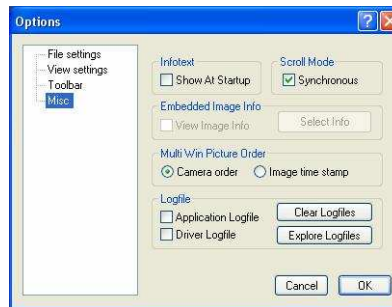

Infotext

Use this command to show or hide the start-up information dialog box.

Synchronous Scrolling

If this item is checked all opened image windows will be scrolled at the same time and in the same manner.

View Embedded Image Info (for pixelfly and sensicam models)

If this item is checked an embedded text will be shown inside the image. See chapter 7.6.

Multi Win Picture Order

With the 'Multi Window' you have the ability to view the images of all active cameras in one window. You can select between 'time' oriented or 'camera number' oriented order. 'Time' oriented means that the sequence of images depends on the individual time stamp of each image. 'Camera number' oriented means that the first image shown is from camera 1, then camera 2 and so on.

Logfile

This option enables the logging of errors and system info. Enable this feature in order to create a support file, which can help to find out troubles with the cameras (see A6 "Customer Service and Trouble Shooting").

Load lookup Table (for monochrome cameras)

With the Pseudo LUT feature you can load any LUT with one of four different formats and you can view the result in the color view window.

Use one of the attached predefined LUTs or define your own.

Direct Record To File


This command is intended for cameras with internal memory. In case of the pco.1300 images are recorded directly onto hard disk by default. With this command you can preset a certain number of images to be stored. If the camera captures images faster than the computer can save to disk, then you will lose images. The displaying of the images doesn't interfere with the record process.

Record with 'Flowfile' (for pixelfly and sensicam models)

Exit

This command exits the program and closes all channel dialog windows.

Window positions, settings and sizes are stored in the windows registry and will be loaded again at next start-up.

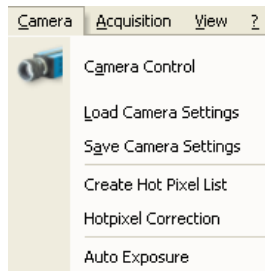

7.4 The Camera Menu

Camera Control

Use this command to open the camera control window (see 6.1).

Load Camera Settings

With this command you can reload camera settings which have been previously saved. You can reload only a file which has been made with the same camera configuration as it was found by the application at start-up.

Save Camera Settings

This command saves the actual camera settings of all existing cameras.

Create Hot Pixel List


Use this command to create a 'Hotpixel-File'. Some pixels tend to become very bright during a long exposure. These pixels are called hot pixel. To remove hot pixels from the image you have to create a hot pixel file first. If you have already a hot pixel file you only have to enable hot pixel correction. The new value of the hot pixel will be calculated from the values of the neighbour pixels. You only can create a hot pixel-file if the last exposure was longer than 5s.

This command will open the hot pixel dialog.

You have to type in the desired count of hot pixels first. Then press the 'Get Hotpixel' button. After this the hot pixel list will be created, using the last acquired image.

Now you have the possibility to save the hot pixel list for later use. All list box items with a checkmark will be saved to the hot pixel file.

You can select or deselect a group of items by highlighting a group and pressing the checkmark in front of one of the highlighted items. Additionally you can open a local menu inside the list box by pressing the right mouse button. You can verify your selection by scrolling through the list and by a comparison to the average value shown in the dialog.

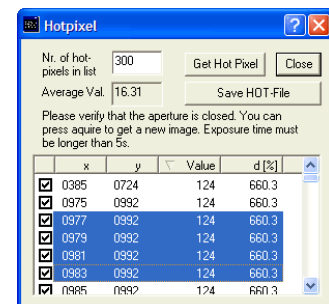

Hotpixel Correction

You can enable or disable the hot pixel correction. If hot pixel correction is disabled, a file dialog will be opened, in which you have to select the correct hot pixel file for the corresponding camera. If you have more than one camera connected you have to select the correct camera number first. If hot pixel correction is enabled this feature will be switched off.

Auto Exposure

Starts some exposures and tries to find an exposure time that captures images with grey values between 1400 and 1500. It re-adjusts the lookup tables in order to display the images with the new settings best.

In case the software is not able to calculate an appropriate setting, this message will be displayed.





7.5 The Acquisition Menu

Live Preview

The 'Live Preview' is useful for fast and easy adjusting and focusing of the camera. The active window will be updated. To see another window, simply click on the window. This option is not available in double shutter mode.

Acquire Picture

Records single images in case of pixelfly and sensicam cameras.

Acquire Sequence

Starts recording images into the system memory according to Trigger Mode selection (see 6.2). During the recording, the camera controls are locked.


Rec. Memory Settings

(for pixelfly and sensicam models)

Recorder Settings

This command opens the Recorder Settings dialog window.

The **Play Settings** are referred to the camera. If more than one camera is connected, the camera number is retrieved from the image window that has received the input focus (is active). If no image window is opened, the dialog is referred to camera 1. If the 'Set to all' checkbox is checked the settings apply to all connected cameras.

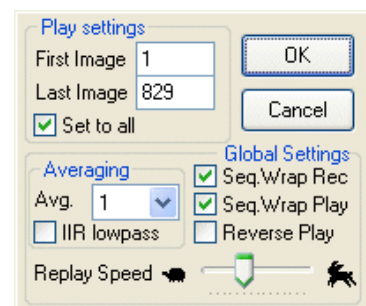

The **Global Settings** are valid for all connected cameras.

Averaging

If in the "Avg." dropdown list a value not equal to 1 is selected, the corresponding number of images is averaged in the buffer, reducing the statistically independent noise.

Another option to reduce the noise is the activation of the "IIR lowpass" IIR-filter: $\text{Image}(\text{act}) = \text{Image}(\text{act} - 1) * 0.5 + \text{Image}(\text{new}) * 0.5$

Replay Speed Slider

The slider can be used to adjust the sequence replay speed, from slow (turtle) to fast (rabbit).

Seq. Wrap Rec

Sequential Wrap Recording sets Recorder Mode (6.6) to "Ring Buffer".

Seq. Wrap Play

The play button starts the display for a sequence as defined by the first image/last image fields and will loop until the stop button is pressed.

Reverse Play

If the "Reverse Play" box is checked, it is possible to play the recorded images in reverse.



7.6 The View Menu


B/W Window

Use this command to open a b/w display window.

Color Window

Use this command to open a color display window.

Convert Control BW

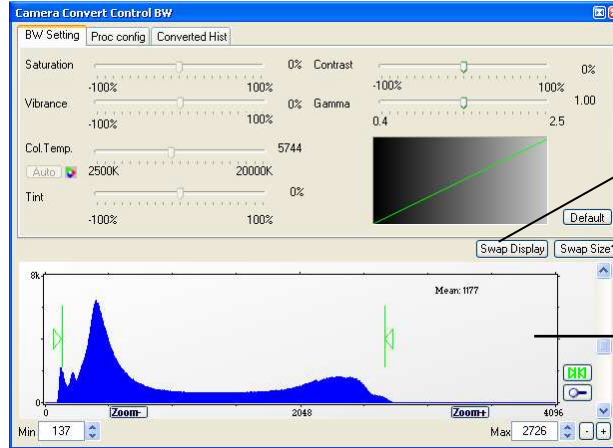


The user can influence how the 12 bit intensity values (x-axis) of the original image are converted into 8 bit values (y-axis) in different ways.

Green sliders in histogram

- left slider = 'Min' controller (corresponds to value "0" of the 8 bit display)
Values below that mark are set to "0", i.e. displayed as black.
- right slider = 'Max' controller (corresponds to value "255")
Values above that mark are set to "255", i.e. displayed as white.

The values in-between are converted into a value between "0" and "255" according to Contrast and Gamma settings. See the small graph, which reflects the calculation.

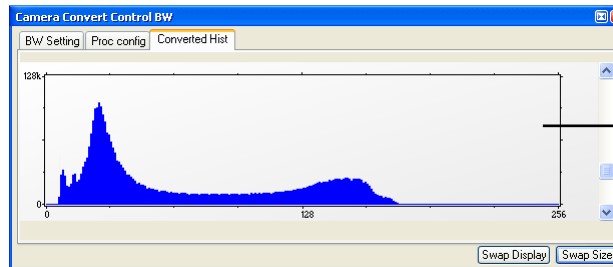

How the Auto Range applies:



Auto Range Peak searches for the minimum and maximum 12 bit intensity values of the image. Given these number the converter scales the 8 bit display (256) within these two values.



Auto Range Crop sets the converter to ignore the extreme intensity values of the image and scales the display in a smaller range. Thus dark or bright light spots, reflections, etc. are cut off.



Min/Max (Automatic) = Continuous Auto Range Crop


7 Camware Features

GPU Processing

In order to benefit from all of Camware's advanced processing features, we highly recommend a NVIDIA CUDA-enabled graphics processing unit (GPU). AMD's ATI graphics processing units are not yet enabled, but will follow soon.

Camware Image Processing

	color	monochrome
Pre Processing		
Noise Reduction	✓	✓
Bayer Processing		
Bayer Demosaicing	✓	
Sharpen	✓	
Refine Process	✓	
Chroma Blurr	✓	
Post Processing		
Offset, Gain (Min/Max)	✓	✓
Color Correction	✓	
Saturation, Vibrance	✓	
Color Temperature, Tint, White Balance	✓	
Gamma, Contrast	✓	✓

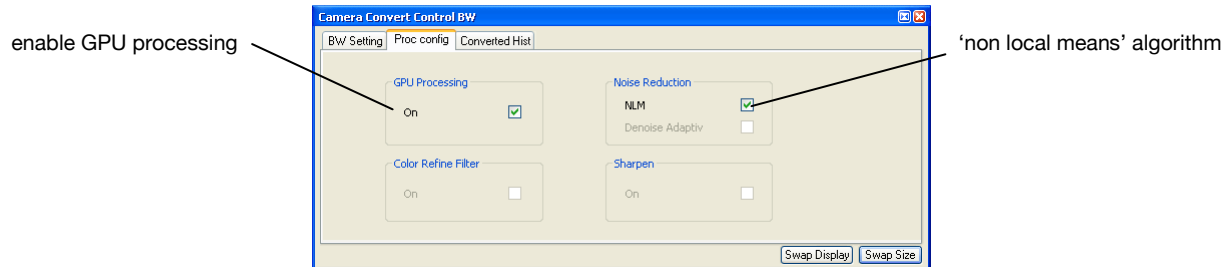


Convert Control Color

This display will be shown only when you connect a color camera to one of your channels.

Browser

The 'Recorder Browser' gives an overview of the images in the Recorder, and helps you in an easy way to select a new current image. Simply click on the desired image to make it the current image. It may then be viewed in the Black/White and Color Window.

The Browser window can show up to 40 pictures per page, thus enabling easy printing of the Browser window. You can switch between the single pages by double clicking inside the Browser window or with the local menu (right mouse button).

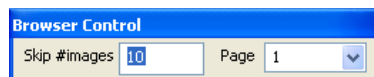


'Skip #images' allows for convenient overview at large numbers of images.

Multi Window

View the images of all active cameras in consecutive order in one window. 'Time' oriented means that the sequence of images depends on the individual time stamp of each image. 'Camera number' oriented means that the first image shown is from camera 1, then camera 2 and so on.

Embedded Image Info (only for pixelfly and sensicam models)

see Timestamp feature in chapter 6.6

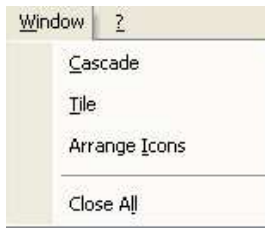

7.7 The Window Menu

Cascade

Show all windows overlapped.

Tile


Tile windows vertically or horizontally.

Arrange Icons

Rearranges all minimized topic icons on the desktop.

Close All

Close all windows.

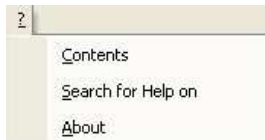

7.8 The Help Menu

Contents

Opens the main page of the online program help.

Search for Help on

Opens an index list for help.

About

This window shows program information and contains helpful support links. In case of a camera problem, support files (see 7.3 Options/Misc/Logfile) may be generated, which can be automatically attached to the email. The support files contain relevant hardware information, the log files and the PCO related registry entries. These files are moved into a zip-archive called "CWSupport.zip", which can be found in the application folder.





7.9 The Local Menu



The local menu is opened by clicking the right mouse button inside the client area of an image window.

Image Properties

Opens an information bubble with main image properties.



Camera Control

Opens the Camera Control window (see chapter 6).

Convert Control BW/Color

Opens the corresponding Convert Control windows (see 7.6).

Zoom

Changes the zoom used to display the image.

Auto Range Peak/Crop

Displays the image in a predefined scaling (see 7.6).

Auto Balance Color

The white balance feature for color cameras.

Auto Exposure

Auto Exposure feature (see 7.4)

Flip

The image will be flipped before display.

Mirror


The selected the image will be mirrored before display.

Rotate

The image will be rotated by 90° clockwise (right) or counterclockwise (left) before it is shown in the image window.

Line Tool

Select 'Line Tool' to show a graph, which shows the intensity of the pixels under the red line shown in the image window. The selection of this menu item determines the location of the first point. After this you can move the mouse to the second point, which can be selected by a 'left click'. Now a 'Line Diagram' opens and the length of the line (unit: pixel) is shown in a bubble and displayed in the window.

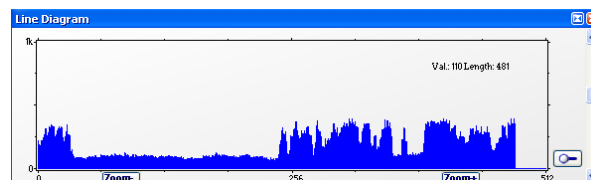

Copy To Clipboard

Copies the active image to clipboard.

7.10 Additional Features

Setting Contrast Area by Mouse

You can control the minimum and maximum values used for the conversion from 12 to 8 bit with the mouse. Move the mouse cursor into a region which should be shown with maximum contrast. Press the shift and the left mouse button. Hold down the mouse button while increasing the size of the rubber band window with mouse moves. After releasing of the mouse button the coordinates of the rubber band window act as a border for calculating the minimum and maximum values.

Setting a new ROI by Mouse

In the same manner you can setup a new region of interest (see ROI in Camera control) for the camera. You only have to press the CTRL (Strg) button instead of the shift button. The coordinates of the rubber band window are used for calculating a new region of interest. You can reset the ROI to maximum by pressing the CTRL (Strg) button and the right mouse button.


White Balance by Mouse

In the same manner you can change the 'white balance'. You only have to press the CTRL (Strg) and the shift button at the same time. The pixel values within the coordinates of the rubber band window are used for calculating a new 'white balance'.

For best results we recommend to use the white balance button in the 'Convert Control Color' (see 7.6) or the 'Auto Balance Color' in the Local Menu (see 7.9) respectively.

Fold Up Window

The Convert Control windows can be minimized/folded up. Just move the pointer over the bar and the window will unfold again.

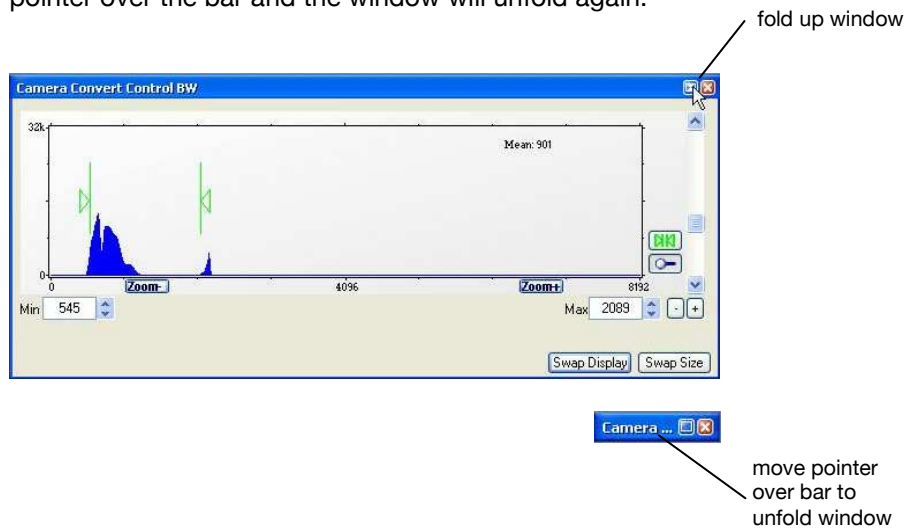

Time Elapse Window

If an image is acquired with an exposure time longer than 5s, the time elapse window will appear to show the remaining time of the current exposure.



Appendix


A1 Technical Data

	unit	setpoint	pco.1300	pco.1300 solar
resolution (hor x ver)	pixel	@extended mode @ normal mode	1424 x 1060 1392 x 1040	
pixel size (hor x ver)	µm ²		6.45 x 6.45	
sensor format	inch		2/3"	
sensor diagonal	mm		11.14	
quantum efficiency (peak)	%	1)	62	13
full well capacity	e ⁻		16 000	
dark current	e ⁻ /pixel · s	@ 10 °C typical	0.05	
image sensor			ICX285AL	
maximum dynamic range	dB	CCD + camera @ 10 MHz	68.5	
dynamic range A/D	bit		12	
read out noise	e ⁻ rms	@ 10 / 20 MHz	6 / 10	
imaging frequency (frame rate)	fps	@ 10 / 20 MHz @ full frame	5.9 / 10.5	5.9 / 11.7
A/D conversion factor	e ⁻ / count		3.5 / 3.8	3.8 / 1.9
spectral range	nm		290 ... 1100	290 ... 1200
exposure time	s		5 µs ... 1 h ²⁾	
anti-blooming factor		@ stand. / @ low light mode (100 ms expos. time)	> 400 / > 4	
smear	%		< 0.002	
optical input			c-mount	
trigger, auxiliary signals		internal / external	software / TTL	
binning (hor x ver)			1x1, 1x2, 2x1, 2x2	
power supply	VDC	typical	18 ... 28 V 24 V	
power consumption	W	Max.	20	
mechanical dimensions (w x h x l)	mm ³		113 x 104 x 110	
weight	kg		1.1	
ambient temperature	°C	range	+10 ... +40	
operating humidity	%	range	10 ... 80	
storage temperature	°C	range	-20 ... +70	
data interface			IEEE1394a	
CE certified			yes	




1) setpoint pco.1300: @ 500nm typical | setpoint pco.1300solar: @ 880nm; +15°C

2) exposure time pco.1300solar @ IR boost: 100ms ... 1h

quantum efficiency

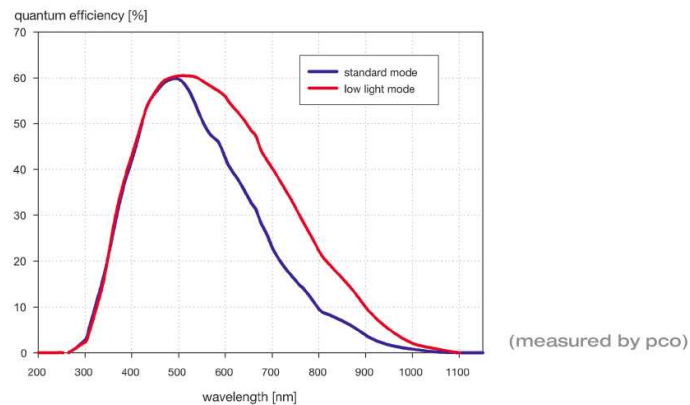

A2 Mechanical Dimensions

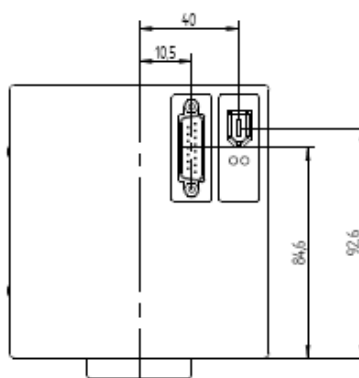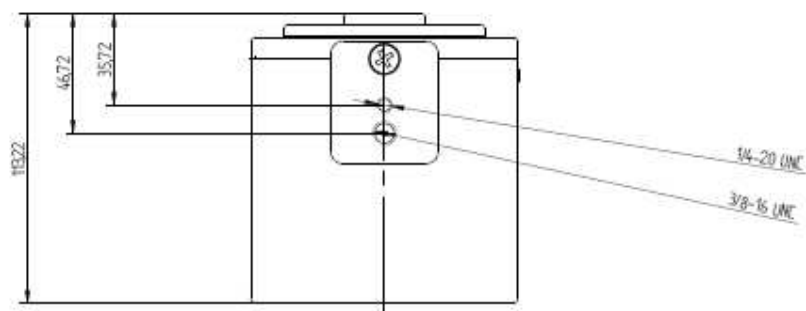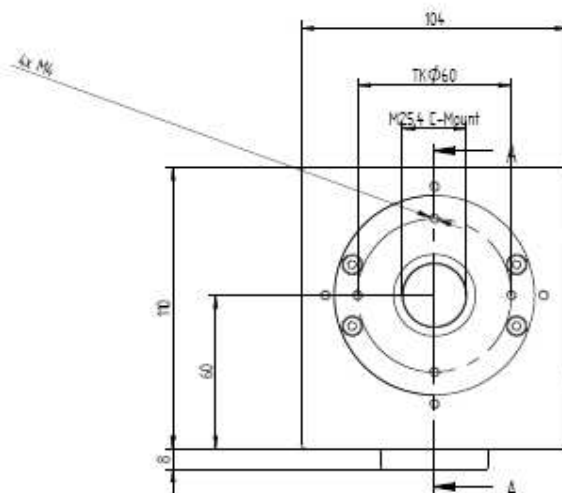

A3 Pinout of 15pin DSUB Connector

PIN	description
1	GND
2	+24V
3	+24V
4	Exp Trig
5	Exp Acquire
6	Busy
7	Exposure
8	GND
9	GND
10	+24V
11	do not connect
12	5VDC/10mA out
13	do not connect
14	do not connect
15	GND

A4 Image File Formats

There are different file formats available for saving camera images with Camware.

b16

The "b16" 16 bit format is similar to the bmp format. However, 16 bit pixel values are used instead of 8 bit pixel values. The 16 bit pixel values represent "unsigned integer values", where the camera (12 bit values) sets the higher 4 bits to zero.

The file format consists either of a "Basic Header" (6 Long-parameter) or of an "Extended Header" (32 Long-parameter), the latter of which is optionally for additional information. There might follow a variable comment field (ASCII code). Finally, there is the actual data set that is saved linearly (as in the case of BMP files).

With the exception of the first value, all parameters are "Long Integers" (4 Byte). The first 6 parameters must always exist. The rest of the parameters, as well as the comment field, are optional.

	Parameter	Function
1	"PCO-"	the first 4 byte are the characters "PCO-"
2	file size	file size in byte
3	header length	header size + comment filed in byte
4	image width	image width in pixel
5	image height	image height in pixel
6	extended header	-1 (true), extended header follows
7	color mode	0 = black/with camera, 1 = color camera
8	b/w min	black/white LUT-setting, minimum value
9	b/w max	black/white LUT-setting, maximum value
10	b/w linlog	black/white LUT-setting, 0 = linear, 1 = logarithmic
11	red min	red LUT-setting, minimum value
12	red max	red LUT-setting, maximum value
13	green min	green LUT-setting, minimum value
14	green max	green LUT-setting, maximum value
15	blue min	blue LUT-setting, minimum value
16	blue max	blue LUT-setting, maximum value
17	color linlog	color LUT-setting, 0 = linear, 1 = logarithmic
18 ... 266	internal use	
Comment file in ASCII characters with variable length of 0..XX. The length of the comment filed must be documented in the "header length" field.		
16 bit pixel data		
	line 1, pixel 1	value of the first pixel
	line 1, pixel 2	value of the second pixel

We recommend that all images should be saved first in the b16 or TIFF format. The advantage is to have the b16 or tiff images available all the time. You will always have the maximum 12 bit information. Please note that not all image analysis programs can accommodate 16 bit data. The 8 bit format saves only the information displayed on the monitor screen. The 12 bit information will be lost and cannot be displayed later.

TIFF

Tag Image File Format, version 6.0 and lower. There is a 16bit monochrome and color image format.

BMP

Windows Bitmap Format, b/w or color 8 bit format-images, which have been saved in BMP format can be loaded later only as 8 bit images, i.e. part of the original information (10bit, 12bit or 14bit) is lost.

FTS

Flexible Image Transport System, Version 3.1. There is a 16 bit image format. The NASA/Science Office of Standards and Technology (NOST) has defined this format. Some programs use the FIT extension for this format.

ASCII

16 bit format, some mathematical programs prefer ASCII data.

JPG

JPEG (named after the Joint Photographic Experts Group who created the standard) is a commonly used method of lossy compression for photographic images. The degree of compression can be adjusted, allowing a selectable tradeoff between storage size and image quality.

JP2

JPEG 2000 is a wavelet-based image compression standard and coding system. It was created by the Joint Photographic Experts Group committee in the year 2000 with the intention of superseding their original discrete cosine transform-based JPEG standard (created 1992).

AVI

Audio Video Interleave is a multimedia container format introduced by Microsoft in November 1992 as part of its Video for Windows technology.

MPG

MPEG-1, similar to JPEG, is a standard for lossy compression of video and audio developed by the Moving Picture Experts Group (MPEG).

WMV

Windows Media Video (WMV) is a compressed video compression format for several proprietary codecs developed by Microsoft. The original video format, known as WMV, was originally designed for Internet streaming applications, as a competitor to RealVideo.

A5 Service and Maintenance

Service

The camera is designed to operate with no need of special adjustments or periodic inspections.



Maintenance

Unplug the camera from any power supply before cleaning it.



Use a soft, dry cloth for cleaning the camera.
Do not use liquid cleaners or sprays.

The lens is best cleaned with pressurized air or with liquid cleaners such as pure alcohol or with special optical cleaners that are available at high quality photo stores.

Never use aggressive cleaning liquids such as gasoline, acetone, spirits or nitro cleanser.

Every time the input window is cleaned, there is the possibility of surface damage. Do not clean the input window unless it is absolutely necessary. Be careful and avoid scratches and damage to the input window surface.

Use a cotton swab dipped in pure alcohol or optical cleaning liquid and wipe only on the glass surface. Do not get any cleaning liquid on the metallic parts such as the lens thread, because tiny detached particles may scratch the surface.

Always store the camera with the protective cap or with a lens mounted to avoid dust and dirt on the input window!



Camera Disposal

The camera includes electronic devices, which can contain materials harmful to the environment. If the camera is to be discarded, please dispose of it in an environmentally responsible manner and use recycling facilities, where available.


A6 Customer Service and Trouble Shooting

If you have a question, which is not adequately addressed in this manual, please contact PCO or your local dealer.

To speed your request, we need the following information:

- Short description of the problem
- Description of your application
- Camera settings
- Type and version of camera software being used
- Camera serial number
- Power supply serial number (pco.power)
- Operating system (PC)
- Processor type (PC)
- Memory
- Graphic card
- Graphic card setup

Please use the email options under the Help menu (see 7.8),

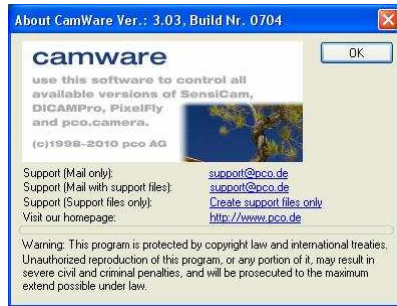

or go to our website: <http://www.pco.de/support-request/>

Before sending the camera for repair, first contact your local dealer or PCO respectively.



When shipping the camera for repair, be certain to carefully pack the camera with proper shipping materials. If possible use the original packaging. Use the protection cap to protect the camera on the lens thread.

A7 Firmware Update

Please contact our support team for further information and assistance!

<mailto:support@pco.de>

<http://www.pco.de/support-request/>



A8 Digital Camera Tools CD-ROM

You find all necessary and some additional software and drivers on the accompanying CD. For the pco.1300 models please select “pco.camera”.



For the latest versions please check our website:


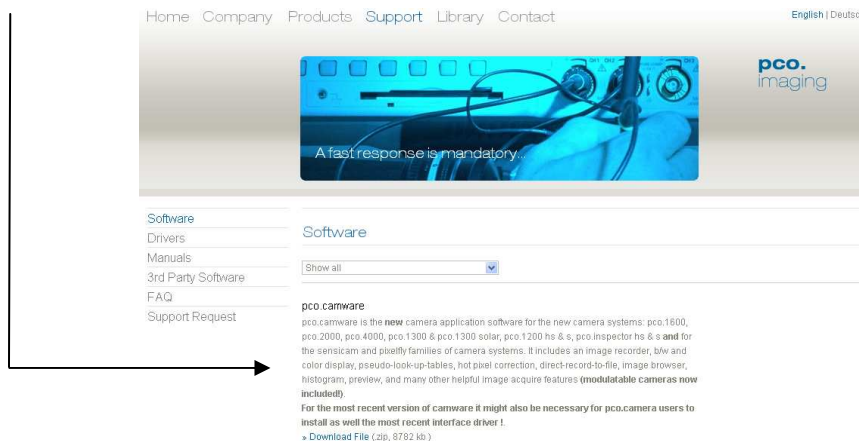


Image Converter

Use the PCO file converter to

- convert 16 bit tif and b16 images to various formats
- display file information
- show thumbnail images

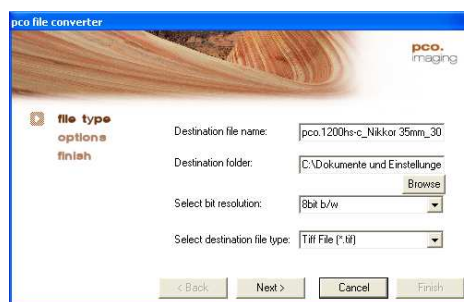

Third Party Image Viewer for PCO b16 images

We recommend the free XnView (www.xnview.com), an efficient multimedia viewer, browser and converter. Please enable “all formats”.

maximize the moment

PCO AG was founded in 1987. The company headquarters in Kelheim employs more than 50 specialists in the development and production of optimized, fast, sensitive camera systems for scientific applications. PCO's range of products includes digital camera systems featuring high dynamics, extremely high sensitivity, high resolution, high speed, and extremely low noise, which are sold in industrial and scientific markets all over the world.

Cameras for every point of view.

The systems produced by PCO AG are cameras and scientific measuring instruments at the same time. Our high-tech systems are mostly the result of manual labor: over 50 highly specialized employees handle development and production at the Kelheim site. We deliver roughly 4.000 cameras a year to customers all over the world. As in every cutting edge technology, dialogue with the user is the main focus of PCO's approach. Worldwide representatives, in cooperation with the in-house marketing division and technical support team, ensure that PCO camera systems are developed in step with the individual requirements of our customers.

pco.
imaging

