


iFLEX® Lasers

High-Performance Laser Systems for Scientific, Analytical and Metrology Applications

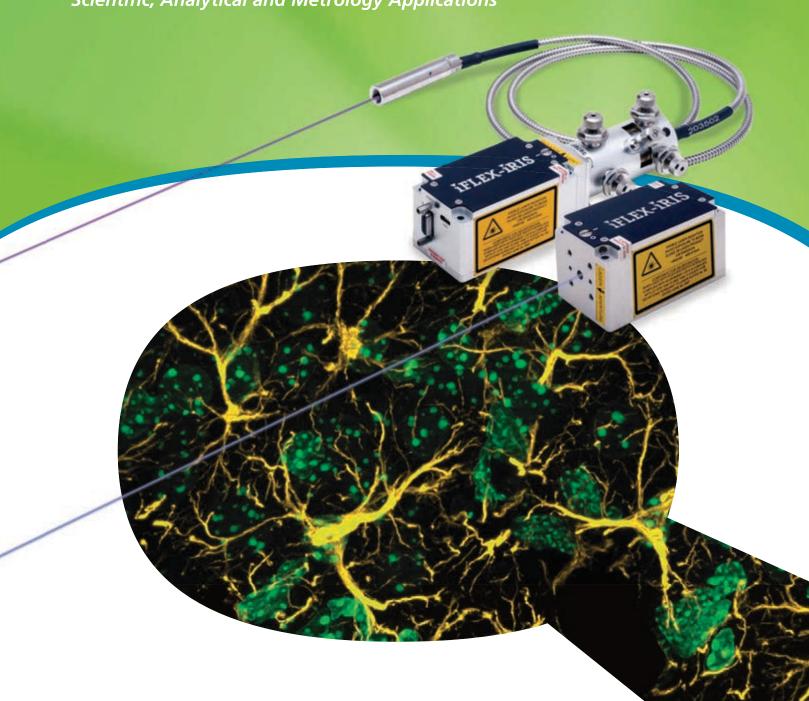

Qioptig, an Excelitas Technologies Company, applications in the areas of medical and life sciences, and aerospace, and research and development.

companies have approximately 5,500 employees in

1877 1898 1966 1969 1984 1991 1996 SPINDLER & HOYER

Rodenstock founded

RODENSTOCK

Spindler & Hoyer founded

Pilkington PE Ltd. founded, which later becomes **THALES Optics**

Gsänger Optoelektronik

founded

Gsänger

Optem International founded

Point Source founded

LINOS

LINOS founded through the merger of Spindler & Hoyer, Steeg & Reuter Präzisionsoptik, Franke Optik and Gsänger Optoelektronik









Content

| iFLEX-iRIS | |
|--|--|
| Introduction & Product Line 04 | |
| iFLEX-iRIS CLM-Concepts and Features 06 | |
| Wavelengths and Specifications 07 | |
| | |
| iFLEX-Gemini | |
| Introduction 08 | |
| iFLEX-Gemini CLM Features 08 | |
| Wavelengths and Specifications 09 | |
| | |
| iFLEX-Viper | |
| Introduction 10 | |
| Wavelengths and Specifications 11 | |
| | |
| Fibers and Optomechanics | |
| Introduction 12-13 | |
| | |
| OEM Expertise & Capabilities 14 | |
| | |
| Total Wavelengths Served 15 | |

2000 2001 2005 2006/2007 2010 2013



Rodenstock Präzisionsoptik acquired by LINOS



AVIMO Group acquired by THALES

Qioptiq founded as THALES sells High Tech Optics Group



Qioptiq acquires LINOS and Point Source as "members of the Qioptiq group"



The new Qioptiq consolidates all group members under one brand



3




Qioptiq is aquired by Excelitas Technologies

Qioptiq designs and manufactures high-performance solid-state laser systems and fiber optics for a range of scientific and industrial applications.

The iFLEX® family of laser technology includes the high performance iFLEX-iRIS™ laser series and the ultra-stable, multi-wavelength laser engine series of iFLEX-Gemini™ and iFLEX-Viper®.

Combine iFLEX lasers with the kineFLEX® single-mode fiber optics manufactured by Qioptiq, to create the world's most stable fiber coupled lasers.

With over 25 years of experience delivering marketleading technology, Qioptiq continues to support customers with demanding applications in semicon, biotech, analytical and industry, through new innovations in iFLEX laser technology.

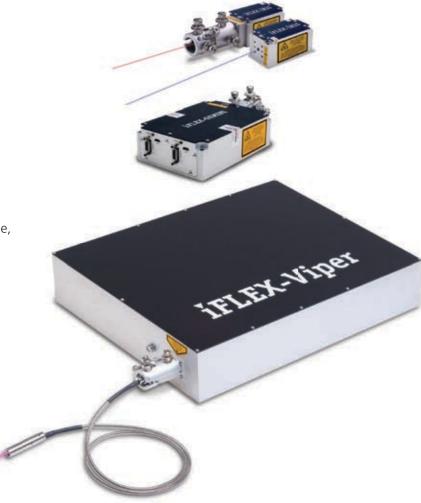

Applications & Features

Features:

- Exceptional power stability
- Unmatched beam pointing stability
- Ultra-low noise performance
- Excellent beam quality
- Fully integrated electronics
- Compact size for easy integration
- End user and OEM systems
- Integrated beam shaping
- "Plug & Play" fiber delivery
- "Set & Forget" alignment

Applications:

- Microscopy
- Flow Cytometry
- DNA Sequencing
- Metrology
- Inspection
- Ophthalmology
- Molecular Imaging
- Dynamic Light Scattering
- Spectroscopy
- Environmental Monitoring



iFLEX-iRIS

Compact, Single-Wavelength Laser Series

The iFLEX-iRIS™ laser series is a range of solid-state, high-performance lasers with low amplitude noise. For ease of use and integration, all wavelengths are offered in the same compact package with the same control inputs. All TEC and smart control electronics are integrated in the laser. They make ideal building blocks for OEM instrument designers and researchers alike.



The innovative Closed-Loop Modulation (CLM) feature allows the lasers to operate with automatic power control feedback in all modes of operation; CW, plus digital, analogue and dual mode modulation. These lasers maintain excellent power stability in all modes of operation and throughout the laser lifetime. Unlike traditional open loop laser modulation, there is no need for laser calibration reset when using iFLEX-iRIS lasers with the CLM feature.

Lasers with CLM are ultra-low noise in terms of RMS, RIN and periodic noise. They also offer precision adjustment at all output power levels. This is very useful for imaging applications where a stable, ultra-low noise source will improve the signal-to-noise ratio and image resolution.



The iFLEX-iRIS lasers can be supplied with a single-mode fiber output. Alternatively, and as requirements change, a single-mode fiber can be added later on by the user to the –X0 laser option, which has the beam exit location centralized in the front face. Thus, the iFLEX-iRIS provides true "Plug and Play" versatility as a free space or fiber

iFLEX-iRIS lasers are designed to fiber couple into the kineFLEX® fiber delivery system. As a result these lasers are ultra-stable when used as free space or fiber coupled. There are standard options for different fiber lengths and either collimated or connector outputs.


A second, free space iFLEX-iRIS laser –X2 option is offered which has the beam exit location offset in the front face, for ease of retrofit in some instruments.

CDRH Compliance for End Users


iFLEX-iRIS lasers are CDRH compliant when used with an iFLEX-iRIS CDRH interlock power supply.





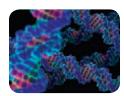

Example: Dual mode modulation

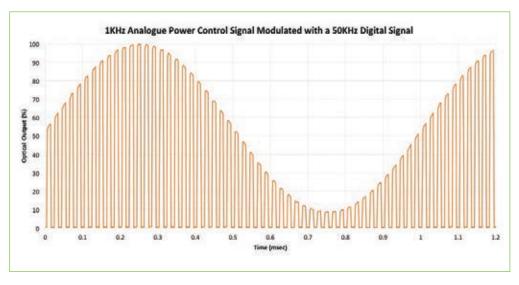

coupled laser.

LASER APERTURE

M3x0.5 IN 4 POSITIONS

iFLEX-iRIS lasers



| | Wavelengths (nm) and Power (mW) | | | | | | | | | | | | | | | | | | | | |
|-----|---------------------------------|-----|-----|-----|-----|-----|-----|-----|-----|-----|-----|-----|-----|-----|-----|-----|-----|-----|----|----|-----|
| 375 | 405 | 413 | 445 | 458 | 473 | 488 | 505 | 515 | 520 | 532 | 561 | 594 | 633 | 637 | 642 | 647 | 660 | 670 | | | 852 |
| 20 | 50 | 100 | 20 | 20 | 75 | 20 | 50 | 20 | 30 | 20 | 20 | 20 | 30 | 20 | 20 | 50 | 80 | 10 | 20 | 70 | 35 |
| 40 | 100 | | 50 | 70 | | 40 | | 50 | | 40 | 40 | | 70 | 100 | 40 | | | | | | |
| 50 | 200 | | 75 | | | 100 | | 60 | | | | | | | 100 | | | | | | |
| 488 | 220 | | | | | 140 | | | | | | | | | | | | | | | |

iFLEX-iRIS fiber coupled lasers

| | Wavelengths (nm) and Power after fiber (mW) | | | | | | | | | | | | | | | | | | | | |
|-----|---|-----|-----|-----|-----|-----|-----|-----|-----|-----|-----|-----|-----|-----|-----|-----|-----|-----|----|----|-----|
| 375 | 405 | 413 | 445 | 458 | 473 | 488 | 505 | 515 | 520 | 532 | 561 | 594 | 633 | 637 | 642 | 647 | 660 | 670 | | | 852 |
| 25 | 30 | 50 | 20 | 45 | 50 | 20 | 30 | 40 | 20 | 25 | 25 | 12 | 20 | 65 | 20 | 30 | 50 | 6 | 10 | 45 | 20 |
| 30 | 50 | | 50 | | | 50 | | | | | | | 45 | | 65 | | | | | | |
| | 100 | | | | | 90 | | | | | | | | | | | | | | | |
| | 130 | | | | | | | | | | | | | | | | | | | | |
| | 150 | | | | | | | | | | | | | | | | | | | | |

iFLEX-iRIS laser specification overview

| Wavelength (nm)** | CLM, 375-520 nm & 633-852 nm | 532 561 594 | | | | | | | |
|--|--|---|--|--|--|--|--|--|--|
| Spatial mode, TEM ₀₀ | $M^2 < 1.2$ | typ | | | | | | | |
| Beam Ø at 1/e ² | 0.7 ± 0.2 r | mm | | | | | | | |
| Pointing stability | < 5 µrad/°C | | | | | | | | |
| Polarization ratio | ≥ 200:1, Vertical ± 2° | | | | | | | | |
| Power supply | 12V DC, | 1A | | | | | | | |
| Base plate temp. | 40 °C maxir | mum | | | | | | | |
| Heat dissipation | 12 W maximum, < | 5W typical | | | | | | | |
| Operation modes | CW / Modulated: Analogue, Digital, Dual Input / Software | CW | | | | | | | |
| Power stability, 8 hrs | < 0.5 % | < 2 % | | | | | | | |
| RMS noise (20Hz - 20MHz) | < 0.05* % | < 0.3* %, <0.1% 561 nm | | | | | | | |
| Peak-peak noise (20Hz to 1MHz) | < 0.5* % | < 3* % | | | | | | | |
| Max periodic noise spike (1KHz -1MHz) | < 0.05* % | <0.3*%, <0.1% 561 nm | | | | | | | |
| CW, power adjustment | 0%, 0.1 - 100% | Off, 50-100%,and 561 nm Off, 15-100% | | | | | | | |
| Digital modulation Bandwidth Extinction ratio Rise / fall time | Digital signal DC to 5 MHz 1,000,000:1 < 100 nsec | OEM options | | | | | | | |
| Analogue modulation Bandwidth Extinction ratio Rise / fall time Power adjustment | 0 - 5V signal DC to 5 MHz 1,000,000:1 < 100 nsec Off and 0.1-100% | OEM options | | | | | | | |
| Dual mode modulation | Two input ports for modulation; same specifications as above. Simultaneous input signals for a) Digital fast On/Off, and b) Analogue power adjustment via external 0-5V input or internal software setting. | OEM options | | | | | | | |
| Communication | micro-USB, RS232 | OEM options | | | | | | | |
| Laser only | 70(L) x 40(W) x 38(H) mm | | | | | | | | |

^{- 20.0} DOWEL SLOT Ø2.02/2.03 x 3 DEEP iFLEX-iRIS Laser FIBER ASSY. 4 MOUNTING HOLES, Ø3.2

iFLEX-iRIS Fiber Coupled Laser

IFLEX-IRIS

4 MOUNTING HOLES, Ø3.2

^{*}Typical performance and wavelength dependant **Center wavelength tolerance typically ± 5 nm.

iFLEX-Gemini Dual-Wavelength Laser Engine Series

The iFLEX-Gemini™ is a series of small, solid-state 2-line laser engines providing a combined, co-axial output beam. It is a turnkey system for OEM instrumentation and researchers. Standard wavelength pairs are listed; custom options available on demand. Applications include: confocal microscopy, flow cytometry, particle sizing, materials testing, optogenetics, metrology, replacement for gas lasers, forensics and medical imaging instrumentation.



Each laser is controlled independently and directly, instead of combining beams through an AOTF. This provides a faster response time as well as instant switching between lines, plus the option for simultaneous emission.

Closed-Loop Modulation Feature

Closed-Loop Modulation (CLM) is offered in the iFLEX-Gemini, for diode wavelengths only. This provides digital




and analogue, and dual input modulation options at DC to 5MHz with high extinction ratios. The CLM feature increases the precision of the power set levels achieved and reduces the noise floor. Thus, it increases the signal-to-noise ratio for many applications.

Fiber Delivery

The iFLEX-Gemini lasers can be supplied with a single-mode fiber output. The user can add or remove the kineFLEX® fiber themselves which provides great flexibility for researchers.

CDRH Compliance for End Users

The iFLEX-Gemini laser engine is CDRH compliant when used with its CDRH interlock power supply option.

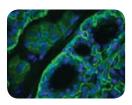

Features & Benefits

Features:

- Output: Combined, coaxial
- Fully independent laser control
- True Off for each wavelength
- Exceptional power stability
- Ultra-low noise performance
- Class-leading beam pointing stability
- OEM and End User options

Benefits:

- No laser alignment required
- Easy to use, portable, turnkey system
- Longer useful lifetime compared to traditional gas lasers
- x10 smaller than equivalent Argon laser
- Direct modulation of each wavelength
- Reliable and repeatable measurements




9

iFLEX-Gemini laser specification overview

| | iFLEX-Gemini | | | | | | | | | | | | |
|-------|---|-----------|----------|----------------|--------|--|--|--|--|--|--|--|--|
| λ1/λ2 | 445 488 515 561 640 647 | | | | | | | | | | | | |
| 405 | 50/50 | 50/50 | - | - | 50/50 | | | | | | | | |
| 445 | 50/50 | | | | | | | | | | | | |
| 488 | 8 50/50 □ 50/30 □ 50/50 | | | | | | | | | | | | |
| | Direct | laser pov | ver (mW) | . Standard λ p | oairs. | | | | | | | | |

| Fiber coupled iFLEX-Gemini | | | | | | | | | | | |
|----------------------------|---|----------|-----------|---------------|-----------|---|--|--|--|--|--|
| (nm) | 445 488 515 561 640 647 | | | | | | | | | | |
| 405 | 30/30 | 30/30 | - | - | 30/30 | | | | | | |
| 445 | - | - | 30/30 | - | - | - | | | | | |
| 488 | 488 30/30 30/20 30/30 | | | | | | | | | | |
| F | iber delive | red powe | r (mW). S | itandard λ pa | irs shown | | | | | | |

| Wavelength (nm) | 405 ± 5 | 445 ± 5 | 488 ± 2 | 515 ± 2 | 561 ± 2 | 640 ± 5 or 647 ± 5 | | | | | |
|---|--|--|---------------------------------------|---|----------------|--|--|--|--|--|--|
| Noise (rms) 20Hz-2MHz | < 0.1* | * % | | < 0.3* % | | < 0.1* % | | | | | |
| Power stability, 8 hrs | | | < | 2 % | | | | | | | |
| Spatial mode, TEM ₀₀ | | M ² < 1.2 typical | | | | | | | | | |
| Laser output beam | | 0.7 mm ± 0.2 mm collimated diameter, collinear | | | | | | | | | |
| Standard fiber options | Type: SM PM fiber Length: 1m, 2m or Output: Collimated | e: SM PM fiber Igth: 1m, 2m or 3m tput: Collimated Ø0.7mm beam or Connector FCP / APC / FCP8 | | | | | | | | | |
| Pointing stability | | < 1 µrad/°C after fiber output < 5 µrad/°C with direct beam (no fiber) | | | | | | | | | |
| Polarization ratio | | ≥ 100:1 | | | | | | | | | |
| Max. base plate temp. | | 40 °C | | | | | | | | | |
| Max. heat dissipation | | | 24 W, < | 5W typical | | | | | | | |
| CW, power adjustment | | 0%, 0.1 | - 100% | | 0%, 15– 100% | 0%, 0.1 - 100% | | | | | |
| Digital modulation Bandwidth Extinction ratio Rise / fall time | | DČ to 1,000 | l signal 5 MHz ,000:1) nsec | | OEM options | Digital signal DC to 5 MHz 1,000,000:1 < 100 nsec | | | | | |
| Analogue modulation Bandwidth Extinction ratio Rise / fall time | | 0 – DC to 1,000 < 100 | OEM options | 0 – 5 V DC to 5 MHz 1,000,000:1 < 100 nsec | | | | | | | |
| Dual mode modulation | Input digital a | N/A | Yes | | | | | | | | |
| Communication | Micro-USB, RS232 OEM option Micro-USB, RS232 | | | | | | | | | | |
| Dimensions | 130 (L) x 90 (W) x 38 (H) mm | | | | | | | | | | |

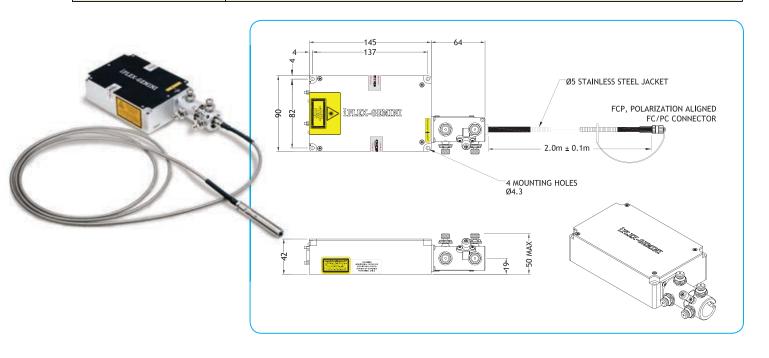

iFLEX-ViperMulti-Wavelength Laser Engine Series

The iFLEX-Viper® is a high-performance, solid-state, multi-wavelength laser engine providing up to 5 lasers in a single system with combined, co-axial output. Robust design eliminates the need for user alignment of the internal laser sources. It is portable and easy to use.

Precision Control

The power adjustment and modulation pattern for each laser is independently controlled; instead of combining beams through an AOTF and allowing the lasers to always emit. Fully independent laser control enables instantaneous switching between wavelengths and simultaneous emission of any wavelength combination. Lasers will only emit when requested, so lifetime may be extended.

Automatic closed-loop control ensures excellent long-term power stability.

The iFLEX-Viper is compatible with a number of commercially available imaging software packages, such as, μ -Manager^{IM} and LabView^{IM}.

Permanent Laser Alignment

Robust, novel, opto-mechanical design in the iFLEX-Viper eliminates the need for user alignment of the internal laser sources. It is a true turnkey system requiring only a drive signal per line to initiate laser emission. The ultra-stable design delivers reliable and repeatable measurements in all applications.



It is a true turnkey system for researchers, easily connected by fiber to microscopes and other instruments. Compact OEM versions also available.

iFLEX-Viper

Fiber Delivery

The iFLEX-Viper is designed to fiber couple into the kineFLEX single-mode, polarization-maintaining fiber delivery system. After the fiber output, the different wavelength beams remain co-axial, polarized and combined as they propagate through any other optics and onto the sample where they overlap. The kineFLEX fiber provides easy connection to microscopes and other analytical and biomedical instruments.

Laser performance is specified and guaranteed after the fiber. The kineFLEX SM PM fibers are offered in different lengths with either collimated or connector outputs.

Features & Benefits

Features:

- Output Beam: Combined, coaxial
- Fully independent laser control
- True Off for each wavelength
- Exceptional power stability
- Class-leading beam pointing stability
- Ultra-low noise performance
- Options: USB, fibers, future upgrades

Benefits:


- Easy to use, portable, turnkey system
- No laser alignment required
- Reliable and repeatable measurements
- OEM or CDRH compliant systems

iFLEX-Viper fiber coupled laser specification overview

| Wavelength (nm) | 640 |) ± 5 | 561 | ± 2 | 532 | 2 ± 2 | 515 | ± 2 | 488 ± 2 | | 445 | ± 5 | 405 | ± 5 |
|--|------------|-----------------------------------|---------|-------|-----------|----------|------------|-----------|-------------|---------|-------------|------------|-----------|----------|
| Nomenclature | I | R | ` | Y | (| 51 | (| Ĵ | | В | | l | ١ | / |
| Power after fiber (mW) | 20 | 50 | 20 | 50 | 20 | 50 | 20 | 40 | 20 | 50 | 20 | 50 | 20 | 50 |
| RYBV - Basic system | | | | | | | | | | | | | | |
| Low power iFLEX-Viper-RYBV | • | | • | | \Q | | \Diamond | | • | | \Diamond | | • | |
| High power iFLEX-Viper-RYBV | | • | | • | | ♦ | | \Q | | • | | \Diamond | | • |
| YGBI - Basic system | | | | | | | | | | | | | | |
| Low power iFLEX-Viper-YGBI | \Diamond | | • | | ◊ | | • | | • | | • | | \Q | |
| High power iFLEX-Viper-YGBI | | \Diamond | | • | | ♦ | | • | | • | | • | | ♦ |
| Noise rms (20Hz – 2 MHz) | | < 0.3 % typ | | | | | | | | | | | | |
| Power stability (8 hours) | | < 2 % | | | | | | | | | | | | |
| Spatial mode, TEM ₀₀ , M ² | | M2 < 1.1 typ, diffraction limited | | | | | | | | | | | | |
| Pointing stability after fiber | | | | | | | < 1 µı | rad/°C | | | | | | |
| Polarization extinction ratio | | | | | | | ≥ 10 | 00:1 | | | | | | |
| Max. base plate temp. | | | | | | | 40 | °C | | | | | | |
| CW power adjustment (per λ) % | 0, 0.1- | -100% | 0, 0.1- | -100% | 0, 0.1 | -100% | 0, 0.1- | -100% | 0, 0.1-100% | | 0, 0.1-100% | | 0, 0.1- | 100% |
| Analogue modulation (per λ) | 0 – | 5 V | 0 – | 5 V | 0 – | 5 V | 0 – | 5 V | 0 – | - 5 V | 0 – | 5 V | 0 – | 5 V |
| Bandwidth | | | | | DC t | o 2MHz, | over 3dE | B bandwi | dth frequ | uency | | | | |
| Dynamic range | | | | | | | ≥ 30 | O dB | | | | | | |
| Rise / fall time over 10 – 90% | | ≤ 350 ns | | | | | | | | | | | | |
| Dimensions laser head | | | | | | 155mm (L | .) x 335m | ım (W) x | 84mm (l | H) | | | | |
| Dimensions controller | | | | 370 | mm (L) x | 322mm (| W) x 85r | mm (H) (c | or H = 91 | mm with | feet) | | | |

[♦] Options for 5th line. Other wavelength combinations are also available.

o Fiber output options: 1m, 2m or 3m lengths, 0.7mm diameter collimated or connectors (FCP, FCP8, APC)

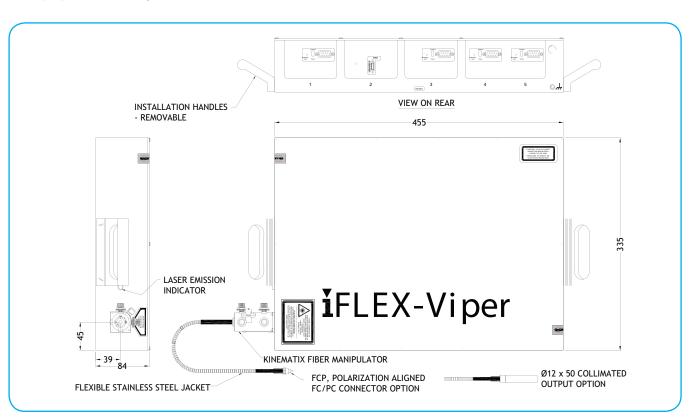

kineFLEXFiber Delivery Systems

The kineFLEX® is a robust SM PM fiber delivery system, suitable for use with most lasers. These single-mode, polarization-maintaining fiber delivery systems deliver the world's best beam pointing stability, making them industry standard in many imaging and precision measurement applications. This fiber delivery system includes: integrated input optics that are pre-focused and optimized for the laser, output optics or connector, and the fiber coupler.

kineFLEX Fibers can be coupled to most lasers:

Single λ : for diode, DPSS, gas etc.

White light λ : 488-640 nm, ArKr, RGB systems

Broadband λ : 400-640 nm, 488-780 nm

ULTRA wide λ : 400-800 nm

High Power: to 500mW standard or custom

Ultraviolet λ: 355 nm



Standard Option Examples

• Wavelengths: 355 nm to 852 nm

Lengths: 1m, 2m, 3m

• Power CW: 100mW or <500mW

• Collimated output: 0.7mm diameter

Connector output: FCP, APC or FCP8



Reasons to use a kineFLEX Fiber System

- Easy beam delivery from "A" to "B"
- Remove hot spots / side lobes / irregularities in the beam profile as the fiber acts as a spatial filter
- World's best pointing stability <1µrad/°C
- Higher power throughput levels
- Fast, and efficient instrument manufacture and servicing with detachable fiber delivery
- Reduce optical errors from multiple interfaces with integrated beam shaping
- Improve instrument stability by removing risk of bulk optic movement through integrating beam shaping optics in fiber
- Robust fiber, safely enclosed laser beam
- Compatible with different environments such as vacuum, UHC, dusty, vibrating
- Custom beam shaping integrated in fiber

OEM Custom Option Examples

- SM, PM, MM & Multi-Channel fibers
- UV-VIS-NIR and lengths to ~40m
- Integrated beam-shaping optics for elliptical, focused, collimated, flat top and parallel spot patterns
- Few Watts into VIS λ SM PM fiber
- Vacuum compatible with options for single core from air to vacuum
- Photonic crystal fiber options

kineMATIX Fiber Coupler

The kineMATIX® is the patented opto-mechanical mount used to align the laser beam into the single-mode fiber. The kineMATIX manipulator is included in the kineFLEX fiber delivery system. A kineMATIX manipulator is included with the kineFLEX Fiber Delivery Systems.

Extremely stable opto-mechanics

It has 4-axes of adjustment (X, Y, tip, tilt) and a centrally located fifth button for easy removal and insertion of the fiber. The design provides sub-micron repeatability and sub-microradian stability. This robust and thermally insensitive design enables the kineMATIX to maintain laser-to-fiber alignment across a wide temperature range and during transit. It truly offers "Set & Forget" laser to fiber alignment.

"Plug and Play" performanace

For over 25 years, the iconic kineMATIX has given our customers confidence to use detachable fiber systems inside OEM instruments and in scientific research. It is the only fiber coupler to offer reliable and repeatable "Plug & Play" performance, since the alignment is retained when the fiber is removed and remains when the fiber is re-inserted.





OEM Expertise and Capabilities

Qioptiq is happy to work with our customers to ensure that each fiber system, laser system or laser engine is optimal for the application, and so we offer custom OEM solutions.


An example of these high-end OEM solutions are the iFLEX-HYDRATM laser engine systems. These are custom designed, multi-wavelength laser engines, with integrated beam-shaping on the fiber output to produce spatially separated beams. These systems effectively combine all the wavelengths needed within the smallest dimensions, and with a SM PM fiber output generating the required beam spot pattern in the flow cell or sample.

They can also be paired with a collection of fiber array to maintain the smallest possible beam paths. This leads to instrument size reduction, with smaller multi-channel laser systems and typically x10 reduction in optics path length.

Engage Flexible Laser Technology™ and true development partnership

With over 25 years supporting demanding applications in industrial manufacturing, biotechnology, clinical diagnostics and semicon, Qioptiq brings design and manufacturing expertise in lasers, fiber optics and multiline laser engines to every partnership.



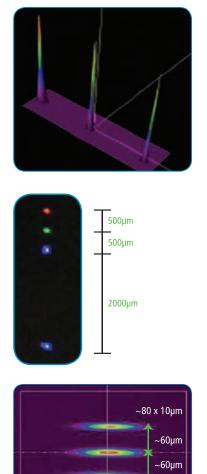

The iFLEX Spectrum

| | Standard Maximum Output Power (mW) | | | | | | | | | | | | |
|-------------------|------------------------------------|--------------------------------------|------------------------------|---|--|--|--|--|--|--|--|--|--|
| | iFLEX-iRIS Laser | iFLEX-iRIS Fiber Coupled Laser | iFLEX-Gemini Laser Engine | iFLEX-Gemini Fiber Coupled Laser Engine | iFLEX-Viper Fiber Coupled Laser Engine | | | | | | | | |
| Wavelength (nm) | | | | | - Carante | | | | | | | | |
| 375 | 50 | 30 | | | | | | | | | | | |
| 405 | 220 | 150 | 50 | 30 | 50 | | | | | | | | |
| 413 | 100 | 60 | | | | | | | | | | | |
| 445 | 75 | 50 | 50 | 30 | 50 | | | | | | | | |
| 458 | 70 | 45 | 50 | 30 | | | | | | | | | |
| 473 | 75 | 50 | | | | | | | | | | | |
| 488 | 140 | 90 | 50 | 30 | 50 | | | | | | | | |
| 505 | 50 | 30 | | | | | | | | | | | |
| 515 | 60 | 40 | 50 | 30 | 40 | | | | | | | | |
| 520 | 30 | 20 | | | | | | | | | | | |
| 532 | 40 | 25 | | | 50 | | | | | | | | |
| 561 | 40 | 25 | 30 | 20 | 50 | | | | | | | | |
| 594 | 20 | 12 | | | | | | | | | | | |
| 633 | 70 | 45 | | | | | | | | | | | |
| 637 | 100 | 65 | | | | | | | | | | | |
| 642 | 100 | 65 | 50 | 30 | 50 | | | | | | | | |
| 647 | 50 | 30 | 50 | 30 | | | | | | | | | |
| 660 | 80 | 50 | | | | | | | | | | | |
| 670 | 10 | 6 | | | | | | | | | | | |
| 730 | 20 | 10 | | | | | | | | | | | |
| 780 | 70 | 45 | | | | | | | | | | | |
| 852 | 35 | 20 | | | | | | | | | | | |
| Measured at laser | beam exit or after | fiber output for fib | er coupled laser sys | items. | | | | | | | | | |



Discover the Q!

Qioptiq delivers cutting-edge technology for all photonic and optical requirements of OEM System development and scientific research alike. Global production capabilities and state-of-the-art manufacturing guarantee an impressive portfolio of products and solutions. Discover the Q for high-performance solid-state laser systems and fiber optics.

Photonics for Innovation

Contact Qioptiq today:

North America +1 (800) 429 0257

United Kingdom +44 (0) 2380 744 500

Europe +49 551 69 35-0

Asia/Pacific +65 64 99 77 77

qioptiq.com qioptiq-shop.com LASERS@excelitas.com

