

Quick Start Guide pco.dimax 3.6 ST series

Required setup

Your computer:

- is equipped with all provided frame grabbers (CLHS interface),
- has the latest version of **pco.camware** (or pco.vision™),
- has the latest version of the correct **PCO Driver** installed.

Your pco.dimax 3.6 ST:

- is powered by the provided power supply and power cable and
- is connected to the computer via all provided FOL cables.

1 The camera switches on once powered up.

2 Start **pco.camware**.

3 Start **Live Preview**.

4 Right-click, check **Continuous Auto Range**.

5 If necessary, adjust the **Exposure time**.

6 Check the image, use your lens or adapter to set the aperture and focus.

7 Once the object is in focus, end **Live Preview** and click **Record**.

