

# Quick Start Guide pco.edge with CLHS interface

## Required setup

### Your computer:

- is equipped with the provided frame grabber (CLHS interface),
- has the latest version of **pco.camware** (or pco.vision™),
- has the latest version of the correct **PCO Driver** installed.

### Your pco.edge CLHS:

- is powered by the provided power supply and power cable and
- is connected to the computer via the provided FOL cable.

**1** Switch on your camera.

**2** Start pco.camware.

**3** Start **Live Preview**.

**4** Right-click, check **Continuous Auto Range**.

**5** If necessary, adjust the **Exposure time**.

**6** Check the image, use your lens or adapter to set the aperture and focus.

**7** Once the object is in focus, end **Live Preview** and click **Record**.

